

# Cell Connect™

The Cell Connect™ is designed to streamline the addition of hematopoietic stem cells and freezing solution into up to four CryoStore™ freezing bags to prepare for cryopreservation.

It is designed to be sterile docked to both your stem cells and freezing media for closed system transfer of products, potentially eliminating the need for a clean room.




## The benefits of using Cell Connect™ include:

- Ease of use saves on training time needed for laboratory personnel
- Efficiency - make 6 connections and you're done
- Closed system reduces contamination issues
- Needle-free processing reduces potential for injury
- PVC tubing allows for sterile docking

# Cell Connect™

## Example of use:

Close all pinch clamps. Attach cells to spike (A) and attached freeze media to spike (B). Attach up to four CryoStore™ freezing bags via male luers (E). Seal off any unused tubes to reduce cell loss. Open pinch clamps to the desired pathway, transfer cells and solutions. A rocker or oscillator is recommended for freeze media addition.

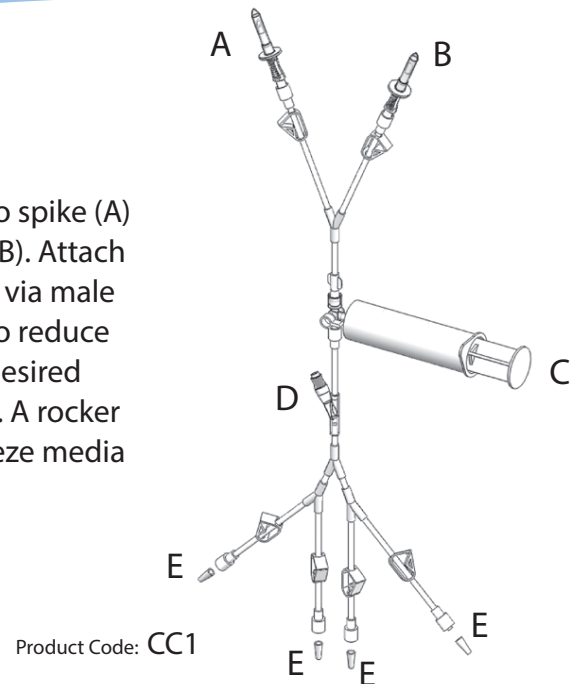

Product Code	Description	Stock/Non-Stock	Case Qty
CC1	Two spikes to a 3-way stopcock, 60ml syringe and needle-free valve to four male luers with 70cm SCD tubing and pinch clamps.	Stock	12
CC2	One spike and male luer to a 3-way stopcock, 60ml syringe and needle-free valve to four male luers with 70cm SCD tubing and pinch clamps.	Stock	12
CC35-5MT	5 male luers with pinch clamps. Intended as an add-on to CC1 or CC2 via sterile dock to create 5 additional terminals for a total of 8 SCD connections for CryoStore™ bag filling.	Non-Stock	12

Individually packed in Tyvek® pouch, supplied sterile, for single use only.

# OriGen

BIOMEDICAL

### Americas & Asia

7000 Burlison Rd, Bldg D  
Austin, TX USA 78744-3202  
Tel: +1 (512) 474 7278  
Tel: Toll Free (800) 233 9014  
Fax: Toll Free (888) 812 8411 or  
+1 (512) 617 1503  
E-mail: sales.us@origen.com

### Europe & Middle East

OriGen Biomedical GmbH  
Romerstrasse 14  
72393 Burladingen, Germany  
Tel: +49-7475-915591  
Fax: +49-7475-914718  
E-mail: sales.europe@origen.com